



Enhancements for best performance with PROGRES GRYPHAX® - Part 2

Additionally to our previously performance enhancement newsletter, please find below enhancements to set-up your customers Laptop or Desktop computer for best performance with PROGRES GRYPHAX® USB 3.0 cameras.

If you have trouble with PROGRES GRYPHAX® USB 3.0 camera connection, please follow the procedure below. You need to deactivate the option: "USB 3 Link Power Management" under advanced power settings from Windows OS (8 / 8.1 / 10) at Power Options.

With the small *.reg file: "Add_USB_3_Link_Power_Mangement.reg" you can display the hidden option: "USB 3 Link Power Management" at Power Options.

Instruction:

1: **Download** the file: "Add_USB_3_Link_Power_Mangement.reg" from link below and run via **double click**.
Note: Administrator permission are necessary to proceed the changes!

https://www.eightforums.com/attachments/add_usb_3_link_power_mangement-reg.47174/

(Note: This is an external link outside our influence area, which can be removed or invalid)


2: Now the hidden option "USB 3 Link Power Management" is visible for all users under category: "USB settings" from advance power settings of "Power Options".

- [-] USB settings
 - [+] USB selective suspend setting
 - [-] USB 3 Link Power Mangement

3: **Open** "Power Options" at "Control Panel" and navigate to option: "Change advanced power settings"
[Change advanced power settings](#)

4: **Deactivate** option: "USB 3 Link Power Management" at advanced power settings. Please see attached screen shot below.

Note: You have to deactivate the options for all power plans separately!

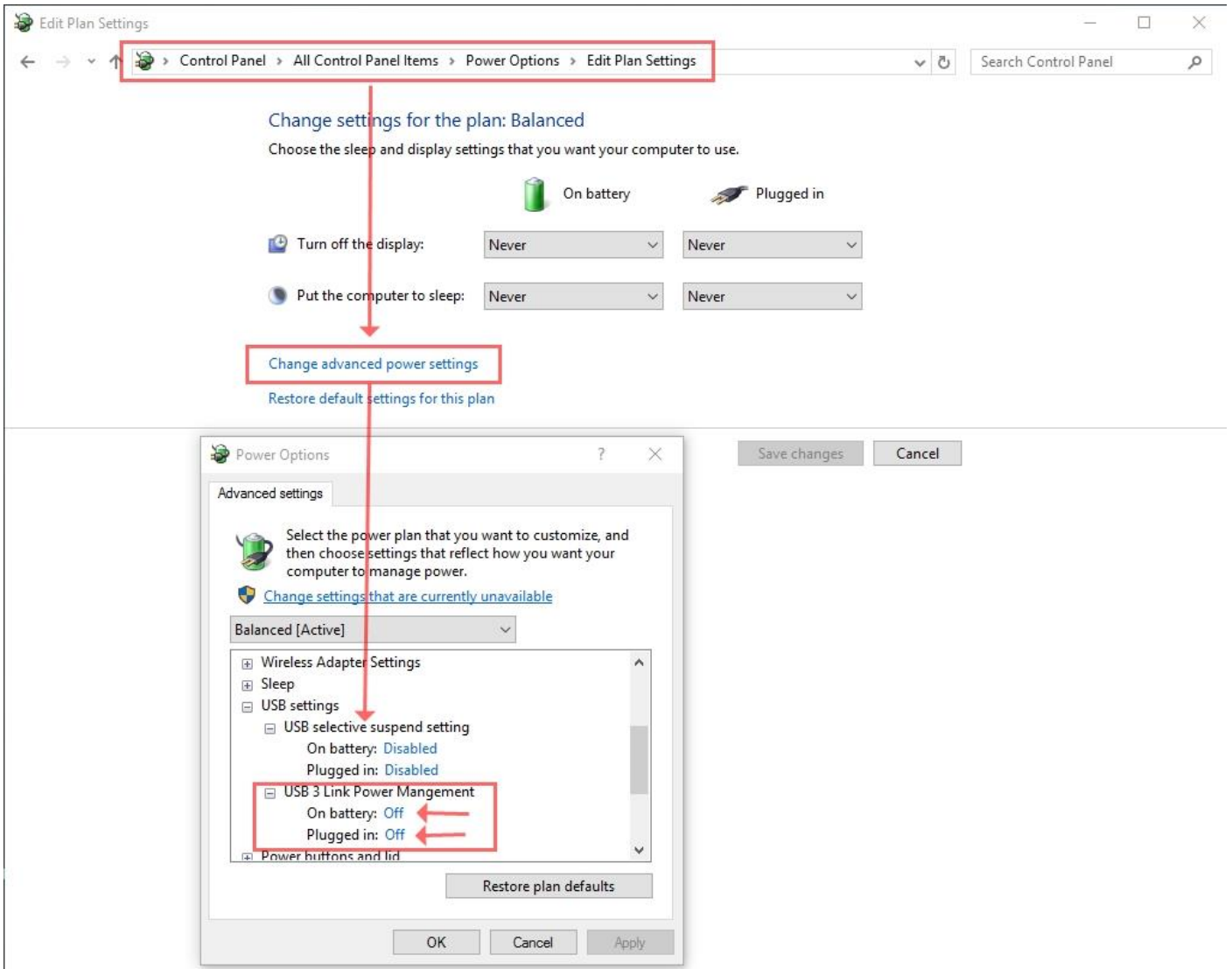

Detailed information about the option: "USB 3 Link Power Management" can be found on the link below:

<http://www.eightforums.com/tutorials/50276-power-options-add-remove-usb-3-link-power-mangement.html>

(Note: This is an external link outside our influence area, which can be removed or invalid)

Limitations:

Option: "USB 3 Link Power Management" not available on Windows 7 or older!

We looking forward to hearing from you and remain sincerely,
Your PROGRES GRYPHAX® Team