




Ease of use –  
giving your work an effortless feel.

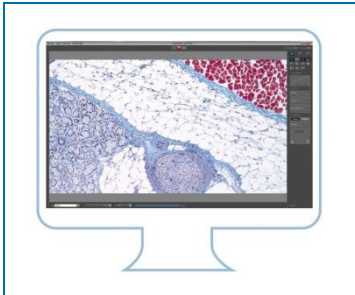

## Shortcuts of PROGRES GRYPHAX® software

PROGRES GRYPHAX® software is workflow-optimized and very easy to use.

Therefore, GRYPHAX software includes a lot of time-saving keyboard shortcuts.

We all want to boost our productivity, so we recommend printing out or saving this list of keyboard shortcuts to enhance your daily work!

Keyboard button:	Shortcuts for:
F1	Open <b>help</b>
F2   ENTER*	Start <b>capturing</b> in "Capture Mode" (single shot, time-lapse, video)
F3	Return to " <b>Live</b> " mode
F5	Activates   deactivates live image <b>freeze</b>
F11	Switch between " <b>Presentation view</b> " and "Standard view"
DEL	<b>Delete</b> selected images from "Gallery" or marked drawings (measurement   annotation)
Z*	Switch <b>zoom</b> between "1:1" mode and "Fit to screen" mode
CTRL** + P	<b>Print</b> selected image(s) from "Gallery"
CTRL** + A	<b>Select all</b> (all thumbnails from "Gallery")
CTRL** + G	Open   close " <b>Gallery</b> "
CTRL** + O*	Open " <b>Storage Options</b> " to change destination folders for gallery (images, videos)
CTRL** + T	Open   close " <b>Toolbar</b> "
CTRL** + D***	Open   close " <b>Treeview</b> " to change and manage media destination folder fast & easy
G	Open   close " <b>Grid</b> " view
ESC*	<b>Leave</b> "Presentation view"   <b>remove</b> drawings during creation (measurement   annotation)
Arrow up   down	Gallery <b>thumbnail</b> up   down
Page up   down	Gallery <b>page</b> up   down
Home	<b>Jump</b> to "Gallery" <b>first</b> page
End	<b>Jump</b> to "Gallery" <b>last</b> page
SPACE BAR	Activates " <b>Magnifier</b> " or " <b>Hand tool</b> "

\*valid as of GRYPHAX version 1.0.14 | \*\*for Mac users: CMD instead | \*\*\*valid as of GRYPHAX version 1.1.6



Learn more about PROGRES GRYPHAX® software and the easy workflows and tools.  
Watch our video tutorials.

[VIDEO TUTORIALS](#)

HELPFUL? We appreciate your feedback.

[FEEDBACK](#)

We looking forward to hearing from you and remain sincerely,

Your PROGRES GRYPHAX® Team

**PROGRES**  
GRYPHAX® SERIES

**JENOPTIK | Healthcare & Industry**  
JENOPTIK Optical Systems GmbH  
Göschwitzer Straße 25  
07745 Jena, Germany